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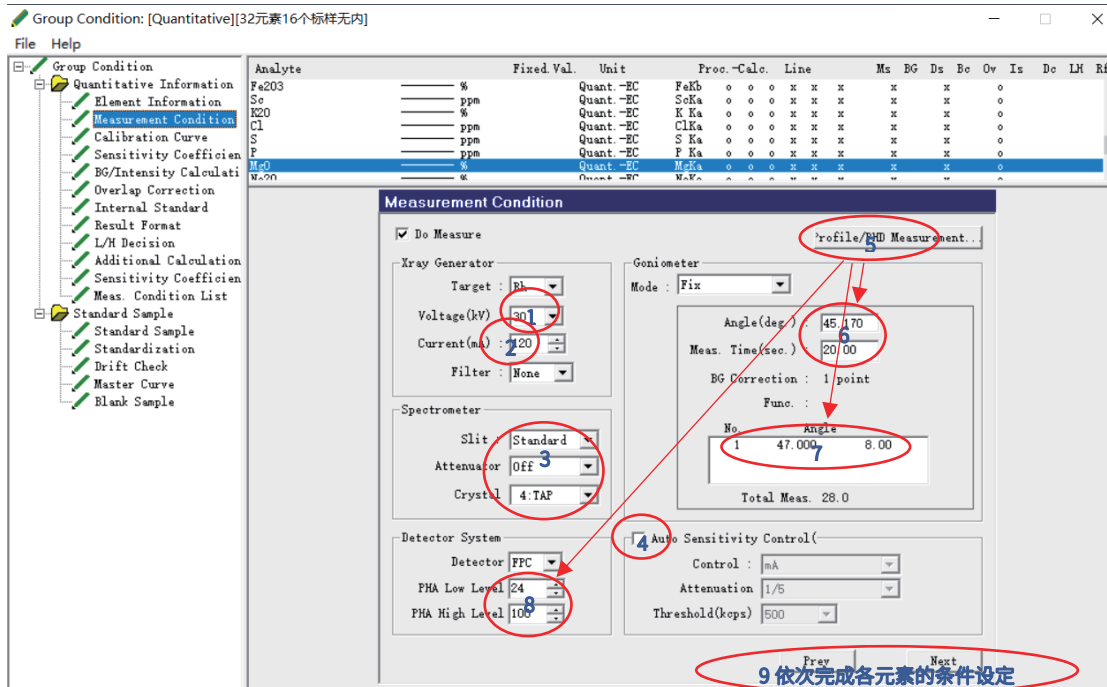
Group Condition: [Quantitative][32元素16个标样无内]

File Help

Group Condition

- Quantitative Information
- Element Information
- Measurement Condition
- Calibration Curve
- Sensitivity Coefficient
- BG/Intensity Calculation
- Overlap Correction
- Internal Standard
- Result Format
- L/H Decision
- Additional Calculation
- Sensitivity Coefficient
- Meas. Condition List
- Standard Sample
- Standard Sample
- Standardization
- Drift Check
- Master Curve
- Blank Sample

Analyte	Fixed Val.	Unit	Proc.	Calc.	Line	Ms	BG	Ds	Bo	Ov	Is	Do	LH	Rf
Fe203	_____	%	Quant.	-EC	FeKb	0	0	0	x	x	x	x	x	0
Sc	_____	ppm	Quant.	-EC	ScKa	0	0	0	x	x	x	x	x	0
K2O	_____	%	Quant.	-EC	K Ka	0	0	0	x	x	x	x	x	0
Cl	_____	ppm	Quant.	-EC	ClKa	0	0	0	x	x	x	x	x	0
S	_____	ppm	Quant.	-EC	S Ka	0	0	0	x	x	x	x	x	0
P	_____	ppm	Quant.	-EC	P Ka	0	0	0	x	x	x	x	x	0
MgO	_____	%	Quant.	-EC	MgKa	0	0	0	x	x	x	x	x	0
MgO	_____	%	Quant.	-EC	MgKa	0	0	0	x	x	x	x	x	0

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Measurement Condition

Do Measure [Profile/PHD Measurement...]

X-ray Generator
Target: Rh
Voltage (kV): 30
Current (mA): 70
Filter: None

Spectrometer
Slit: Standard
Attenuator: Off
Crystal: 1-LiF

Detector System
Detector: PFC
PHA Low Level: 20
PHA High Level: 80

Profile/PHD
File View

Group Condition
Aperture: 30.0
Atmosphere: Vacuum
 Spin
Rotation: 0
 Low Speed Evac./Purge
 Vacuum Stabiliz.
 Higher Order Prof.: Pressure to 30

Channel Condition
Target: Rh
Voltage: 30
Current: 70
Filter: None
Slit: Standard
Attenuator: None
Crystal: 1-LiF
Detector: PFC
PHA Low Level: 20
PHA High Level: 80

Measurement
Position: 1
Measurement: 2theta
Start Stop
Eject

2theta Condition
Scanning: Fixed Angle Step
Peak Angle: 97.75 Step Angle: 0.050
Start: 65 Speed: 8.0
End Angle: 101 Time: 0.375

PHD Condition
Peak Angle: 97.75
Start: 0 Step Width: 2
End: 200 Time: 1

Profile Monitor: [2theta]

View Process Data Set Measurement Angle

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Angle = 62.520 deg. Intensity = 14.968

Measurement

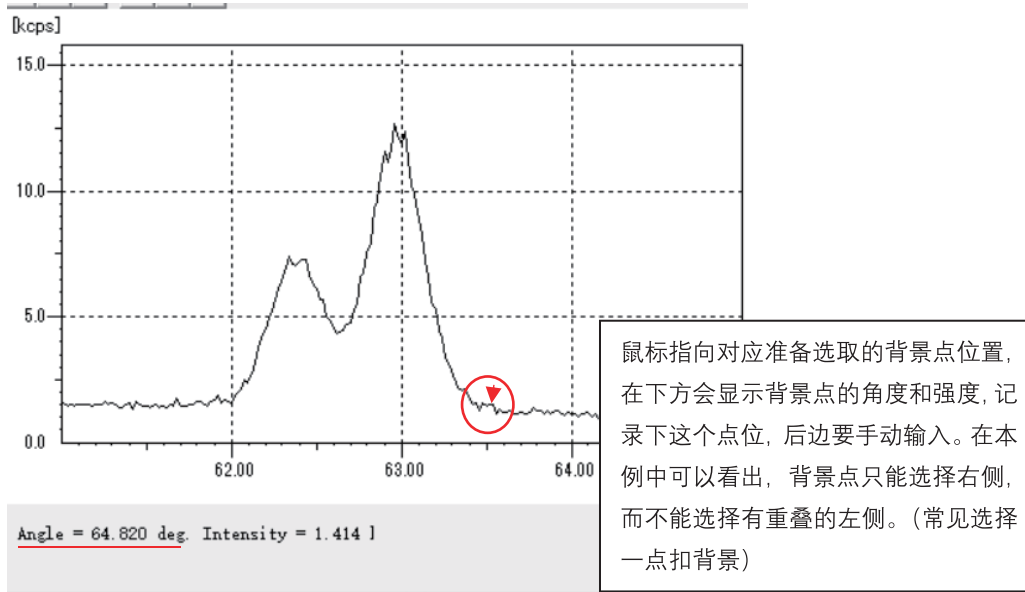
Position: 8
Measurement: 2theta PHD Intensit
Start Stop
Eject

2theta Condition
Scanning: Fixed Angle Step Continuous
Peak Angle: 62.97 Step Angle: 0.020
Start: 61 Speed: 8.0
End Angle: 65 Time: 0.15

PHD Condition
Peak Angle: 62.97
Start: 0 Step Width: 2
End: 200 Time: 1

OK

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Group Condition: [Quantitative][32元素16个标样无内]

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